

GP-105 pH MEASURING HEAD FOR pH METERS OF 100 SERIES

GP-105 may be used with all meters with pH measuring function from the 100 series. Those are: CP-101, CP-102, CP-103, CP-105 and CX-105.

The GP-105 head is used for pH and temperature measurements of samples without or with only small amount of deposits. The head can't be used for viscous samples or those with deposits as they may clog the diaphragm.

The head may be used for measurements of:

- drinking water,
- tap water,
- samples of water from rivers and lakes,
- swimming pool water
- rain water etc.

In the plastic housing there are separate glass measuring and reference electrodes and temperature sensor.

Replacing the head is trouble-free, it is necessary to fit special shape of the head to the meter's body and tightly screw the nut.

The electrodes should be kept active by keeping a sponge, placed in the protective cap, wet. For this purpose distilled water should be used.



TECHNICAL DATA

Measuring pH range	0 ÷ 14.00 pH
Measuring temperature range	-5 ÷ 60 °C
Working temperature range	-5 ÷ 60 °C
Outer housing	ABS
Diaphragm	ceramic
Membrane	glass

ELMETRON

41-814 ZABRZE ul. W. Witosa 10 POLAND
tel. +48 32 273 81 06, fax +48 32 273 81 14

info@elmetron.com.pl; www.elmetron.pl